




No front panel should be a hostage!

Version: 0.61.3

Getting Started Guide:

## Downloading WebPager

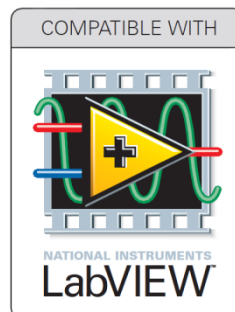


January 2015


Let's see how to download the WebPager software. It will be done in five steps. If in doubt, watch [video tutorial 1](#) on [documentation page](#).


**Step 1:** [Download 'WebPager Downloader'](#) first. Download link is presented on the homepage:



After pressing 'download' your browser will start downloading an OGP installer with WebPager downloader (for example,  [vfp\\_soft\\_com\\_webpager-0.61.3-0.ogp](#)).

**Step 2:** Launch VI Package Manager (VIPM) from JKI. If you don't have VIPM installed on your computer, you can [download it for free from JKI website](#). In VIPM you should choose LabVIEW version and bitness. In this example we choose LabVIEW 2012, 32 bit:

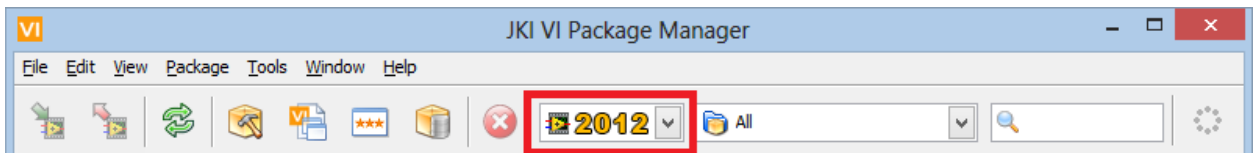

**Step 3:** Run (with VIPM) the OGP installer you downloaded in step 1 (e.g.,  `vfp_soft_com_webpager-0.61.3-0.ogp`).

This is one of the two standard LabVIEW package formats that can be opened with VI Package Manager (VIPM) from JKI. If you don't have VIPM installed on your computer, you can [download it for free from JKI website](#).

**Step 4:** Install WebPager. You can read about this in the next [getting started tutorial 'Installing WebPager'](#).

Don't forget to go through other documentation if you need it.

<https://vfpssoft.com/webpager/documentation>

We are improving this product on a daily basis and value any feedback you could provide. So, if you have any questions, concerns, bug reports, feature requests, propositions, etc. – please drop us a line to [webpager@vfpssoft.com](mailto:webpager@vfpssoft.com).

**LabVIEW**ishly yours,

**WebPager team.**

No front panel should be a hostage!